Apéndice técnico
Technical Information

Volumetría

Volumetry

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Accuracy classification

Volumetric glassware are divided into two groups (classes): class A and class B. The class of the volumetric glassware indicates the degree of accuracy and suitability for official certification. And also the delivery and waiting times. All tolerances fulfil the limits of UNE and ISO standards.

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Delivery and waiting times

The delivery time is the time needed for the meniscus to pass from the upper mark to standstill which is mostly the tip of the pipette or burette.

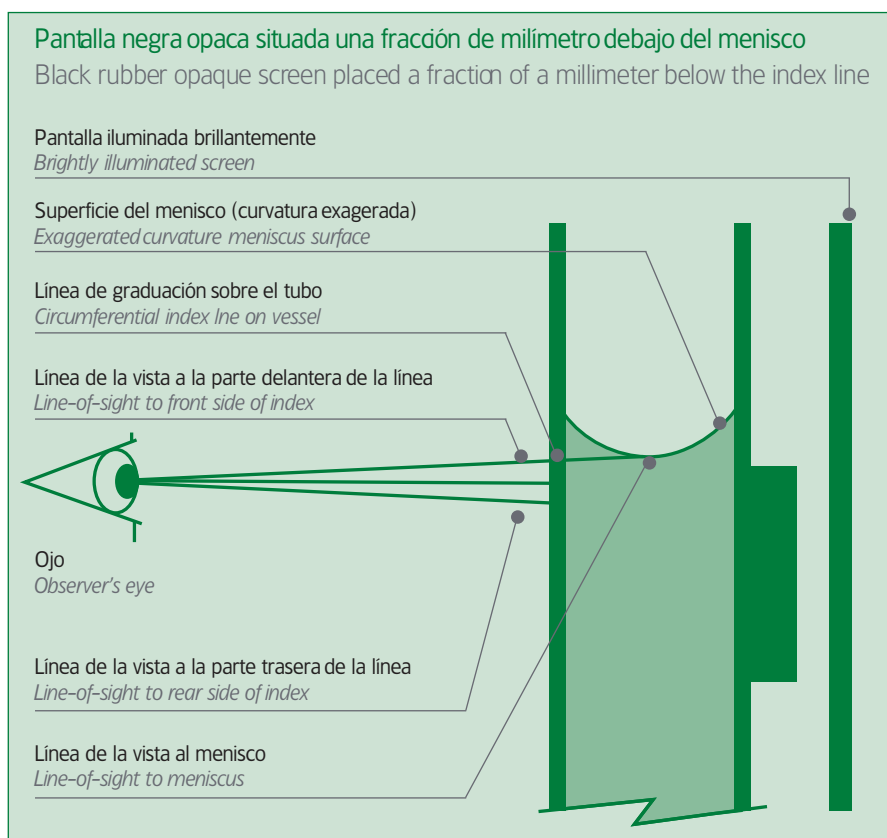
The waiting time starts after the delivery process finishes. During the waiting time remaining liquid continues to flow down from the glass wall. For the most accurate measure of volume with burettes and pipettes someone has to pay attention to the waiting time. Specially with pipettes you have to wipe off the pipette tip on the wall of the glassware that is used.

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Measuring the volume

ISO Standard requires the volume to be read at the lowest point of the liquid level. This lowest point has to touch the upper limit of the graduation mark. Meniscus is the curvature in the surface of the liquid. The curvature is caused by the surface tension of the liquid. On the volumetric glassware the shape of the meniscus is concave (curved downwards). For accurate and most precise results the eye of the observer must be at the same level of the meniscus. To get a better, faster and precise reading and measurement a black strip should be held at the back, just below the graduation mark and meniscus line.

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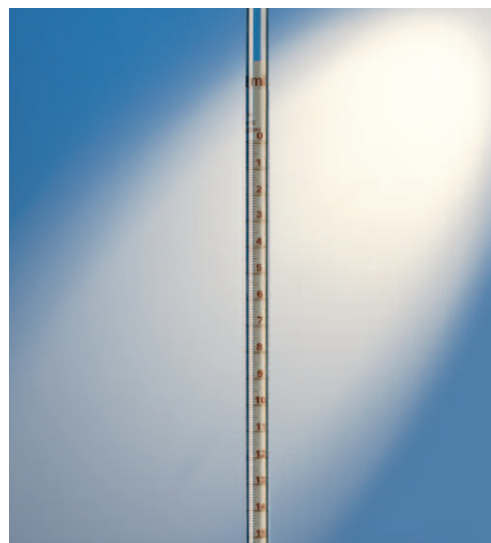
Schellbach stripe

The Schellbach stripe is a narrow blue line on white stripe at the background. Schellbach stripe is printed at the back of some volumetric products such as burettes for easier reading. The light fraction makes two arrows head points to appear at the meniscus line. And the reading point is exactly where these two arrows head meet each other.

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How to use volumetric flasks

Volumetric flasks are mainly used for the preparation of highly accurate standard solutions. They are calibrated according to "to contain". First put the exact and precise weight of the substance into the flask. Then fill the volumetric flask with water (distilled water) about half of the volume and mix the content fully. Then fill in with more liquid just a bit below of the calibration ring mark. Then add some more water so that the meniscus will be exactly at the same level of the ring mark. Pay attention to read the meniscus at the eye level and pay attention not to wet the neck wall above the ring mark level.

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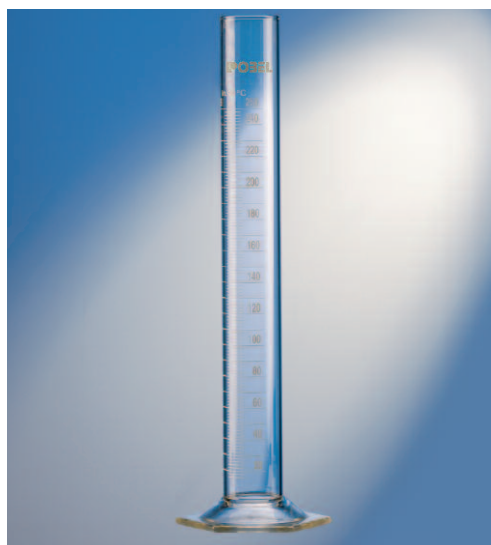
How to use cylinders

Cylinders are mainly used for exact measurement of liquids and they are calibrated "to contain". Fill in with liquid so that meniscus will be exactly at the same level of the desired mark. Pay attention to read the meniscus at the eye level.

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How to use burettes

Burettes are used for titration and are calibrated "to deliver". First rinse the burette very well with the solution you will use for titration. Then fill the burette a bit higher than the zero mark. To activate the stopcock drain the solution but never below the capacity of zero mark. Pay attention that there will be no air bubbles left inside the burette. Fill more titration solution up to max 10mm above the zero mark and pay attention not to wet the burette wall above the liquid level. By reading the meniscus at your eye level drain the solution to the exact level of the zero. At this step you have to wipe the drops at the tip of the burette tip. Then you have to open the stopcock very slowly and start the titration and also pay attention not to touch the tip to the wall of the container you use. When the colour changes stop the titration process and wait for 30 seconds for the waiting time. You have to read the discharged volume at your eye level. Do not forget to wipe the remaining drop of liquid at the burette tip since this is also a part of the titrated volume.

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How to use pipettes

Pipettes are calibrated "to deliver". The first stage is to fill the pipette: for filling process you always have to use a pipette filler or a pipette pump. Fill the liquid max 10mm above the volume mark. Then you must wipe and make dry the outer surface of the pipette. Hold the pipette to your eye level and place the tip of the pipette to the wall of a vessel with a slight angle. Then you have to discharge the excess volume until the meniscus reaches the calibration mark of the pipette. Second step is to empty the pipette. You must empty the pipette by placing the tip of the pipette to the wall of the vessel with a slight angle. Pay attention to empty the pipette completely (or until the 2nd mark of bottom graduation line).

Propiedades del vidrio

Properties of glass for laboratory

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The manufacture of the glassware must reach some physical and chemical properties that make it suitable for use. The one which has been consolidated over the years as a suitable glass is the "borosilicate glass 3.3."

Physical properties

The use in the lab requires a satisfactory product in the range of temperature and pressure. The borosilicate glass POBEL, due to its high content of silica it has been possible to achieve the lowest linear coefficient of expansion ($3.3 \cdot 10^{-6}$). Thanks to this property glass remains indifferent to sudden changes in temperature as well as to thermal shock, depending, of course, on the thickness of the walls and the shape of the product.

However, it is not advisable to heat the glass pieces above the 450 °C so that a permanent tension can occur in the glass. In case that the temperature exceeds 450 °C during handling it is desirable to leave it to cool for several hours at room temperature.

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Chemical Properties

The borosilicate glass POBEL does not contain heavy metals or zinc or arsenic and further its chemical indifference shown in the fact that after repeated autoclave sterilization drastic exceeding 150 °C does not happen any damage on the surface of the glass.

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The pieces of borosilicate glass POBEL are very resistant against the influence of water and acids except the hydrofluoric acid and phosphoric acid concentrated acids while they are in use, because at the same time form a layer on the surface of silica.

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PTFE is the most important member of this family. It has highly stability against most chemicals and it is resistant even over 260 degrees Celsius. Another members of this family are PFA and FEP.

Very high melt viscosity of PTFE prevents the use of usual processing techniques for thermoplastics. Like in powder metallurgy, PTFE material is first preformed usually at room temperature, then sintered above the melting point (typically 370°C), and then cooled. The thermal degradation is negligible in the processing. Additives are seldom used with PTFE.

The molecular configuration of PTFE brings to its surfaces a high anti-adhesiveness.

PTFE is one of the most thermally stable plastic material. There are no appreciable decompositions at 260°C, so that PTFE, at this temperature, still possesses the greater part of its properties. Appreciable decomposition begins at over 400°C.

TetraFluorEthylene-Perfluorpropylene(FEP) is produced by copolymerization of tetrafluoroethylene and hexafluoropropylene. It is a relatively soft thermoplastic with lower tensile strength, wear resistance, and creep resistance than many other engineering plastics. However, it is chemically inert and has a low dielectric constant over a wide frequency range. FEP possesses a very high degree of stress crack resistance, a low coefficient of friction, exceptional dielectric properties, heat resistance, retention of properties after service at 204°C with useful properties at -270°C, and meets FDA 21CFR.177.1550.

FEP has high transparency (with good transmittance of UltraViolet and visible wavelengths.) It has long term weatherability and excellent resistance to ozone, sunlight and weather. FEP offers the lowest refractive index of all thermoplastics with low light reflection (the same as water).

PerFluoroAlkoxy (PFA) offers similar properties to FEP, but is considered more of a premium resin. PFA is preferred when extended service is required in hostile environments involving chemical, thermal, and mechanical stress. PFA offers high melt strength, stability at high processing temperatures, excellent crack and stress resistance, a low coefficient of friction, and more than 10 times the Flex life of FEP.

It has high resistance to creep and retention of properties after service at 260°C, with useful properties at 95°C. PFA also meets FDA 21CFR.177.1550.

CARACTERÍSTICAS FEATURES	PTFE	PFA	FEP
Aspecto / Appearance	Opaco / Opaque	Traslúcido / Translucent	Claro / Clear
M.p. °C	327	305	270
Máx. temp. uso continuo °C / Max. temp. continuous use °C	260 - 290	250	205
Máx. temp. uso intermitente °C / Max. temp. intermittent use °C	315	280	230

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Before sterilizing any items of plastic laboratory ware, verify that no contamination or residues are present. Their presence could destroy plastics during sterilization or autoclaving. Observe the tolerated temperature ranges of plastics when autoclaving plastic laboratory ware. Remove any stoppers, fittings, or caps from plastic laboratory ware prior to autoclaving. Plastic vessels should be autoclaved separately from their closures and other fittings. Autoclaving vessels with their closures in place can lead to deformation and destruction of the vessels.

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This data has been collected from chemical resistance tables supplied by manufacturers and as such serves as guideline for chemical resistance. The actual resistance towards a certain chemical depends always on conditions applied.

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DIN abbreviation, heat and chemical resistances of the plastics

ABR. DIN ABBREVIATION	DENOMINACIÓN POLYMER NAME	TEMPERATURA DE USO TEMPERATURE OF USE		MICRO-ONDA* MW*	AUTOCLAVABLE AUTOCLAVE
		DESDE FROM	HASTA UP TO		
ABS	Copolímero de acronitrilo - butadieno - estireno Acrylonitrile-butadiene-styrene copolymer	-40°C	+85 (100)°C	✓	✗
E - CTFE	Etileno - Clorotrifluoroetileno Ethylene chlorotrifluoroethylene	-76°C	+150(170)°C	✓	✓
ETFE	Etileno - Tetrafluoroetileno Ethylene tetrafluoroethylene	-100°C	+150(180)°C	✓	✓
FEP	Tetrafluoroetileno - Perfluoropropileno Fluorinated ethylene propylene	-200°C	+205°C	✓	✓
HDPE	Polietileno de alta densidad High density polyethylene	-50°C	+80 (120)°C	✓	✗
LDPE	Polietileno de baja densidad Low density polyethylene	-50°C	+75 (90)°C	✓	✗
MF	Melamina Melamine formaldehyde		+80 (120)°C	✓	✗
PA	Poliamida (PA6) Polyamide	-30°C	+80 (140)°C	✓	✗
PC	Policarbonato Polycarbonate	-100°C	+135(140)°C	✓	✓
PE	Polietileno (véase HDPE / LDPE) Polyethylene (see HDPE/LDPE)				
PFA	Perfluoroalcoxy Perfluoroalkoxy	-200°C	+260°C	✓	✓
PMMA	Polimetilo metacrilato Polymethylmethacrylate (acrylic)	-40°C	+85 (90)°C	✓	✗
PMP (TPX®)	Polimetilo penteno Polymethylpentene	0°C	+120(180)°C	✓	✓
POM	Poliosimetileno Polyoxymethylene	-40°C	+90 (110)°C	✗	✗
PP	Polipropileno Polypropylene	-10°C	+120(140)°C	✓	✓
PS	Poliestireno Polystyrene	-10°C	+70 (80)°C	✓	✗
PTFE	Politetrafluoroetileno Polytetrafluoroethylene	-200°C	+260°C	✓	✓
PVC	Polivinilcloruro Polyvinyl chloride	-20°C	+80°C	✓	✗
PVDF	Fluoruro de polivinilo Polyvinylidene fluoride	-40°C	+105 (150)°C	✓	✓
SAN	Estirol - acrilonitrilo Styrene acrylonitrile	-20°C	+85 (95)°C	✓	✗
SI / NR	Caucho de silicona / Caucho natural Silicone rubber / Natural rubber	-50°C	+180(250)°C	✓	✓

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* MW - Micro-wave. Take note of the temperature resistance.

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General table of resistance of polymers against different substances

GRUPOS DE SUSTANCIAS A +20°C SUBSTANCES AT +20 °C	Plástico						PTFE PFA FEP			
	LPDE	HDPE	PP	PMP TPX®	PS	SAN		ECTFE ETFE	PC	PA
Alcoholes, alifáticos Alcohols, aliphatic	A	A	A	A	A	A	A	A	B	L
Aldehidos Aldehydes	B	B	B	B	N	N	A	A	L	L
Bases Bases	A	A	A	A	A	A	A	A	N	L
Ésteres Esters	B	B	B	B	N	N	A	A	N	A
Hidrocarburos, alifáticos Hydrocarbons, aliphatic	L	B	B	L	N	N	A	A	L	A
Hidrocarburos, aromáticos Hydrocarbons, aromatic	L	B	L	L	N	N	A	A	N	A
Hidrocarburos, halógenos Hydrocarbons, halogen	N	L	L	N	N	N	A	A	N	B
Cetonas Ketones	B	B	B	L	N	N	A	B	N	A
Agente oxidante (ácidos) de alta intensidad Oxidizing agent (acids), high intensity	L	L	L	L	L	N	A	A	N	N
Ácidos, diluidos, de baja intensidad Acids, diluid, low intensity	A	A	A	A	B	B	A	A	A	L
Ácidos, concentrados, de alta intensidad Acids, conc., high intensity	A	A	A	A	B	L	A	B	N	N

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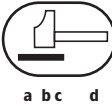
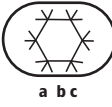
The gloves for specific uses are according the standards detailed below.

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“ CE - Only for low risks ”

It is clearly defined as low risks the gardening, for example, or temperatures from 40 up to 50 °C.
The expression “ CE - Only for low risks ” is marked on the glove

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1 = Very good
2 = Good
3 = Acceptable
4 = Poor

This table provides only general information. The resistance of a glove is influenced by factors such as temperature, concentration of chemicals, thickness of the glove and soak time. To be used is recommended a special pre-test.